

Environmental Shaker Incubators

Dual Action Shaker, MaxQ 2508 (PSC 555, NSC 338, 340, 448)

Every time you use an Incubator Shaker, you MUST sign the logbook.

Use only short screws.

Secure all tubes and flasks – they shouldn't move when shaking.

If you set the temperature for anything besides 37°C, please leave a note and change it back when you are done.

If you hear any kind of grinding sound, STOP.

Contact: Sonja Young (404) 413-5363; sstovall@gsu.edu

MaxQ 2508 Dual Action Shaker (PSC 555, NSC 338, 340, 448)

1. Turn shaker on – front panel in the center.
2. Turn speed control clockwise to increase speed and counterclockwise to decrease speed.
3. Timed shaking. Rotate timer past the 5-minute setting and then adjust the shaking speed. After shaking motion is started, rotate the timer to any timed interval from 1 minute to 60 minutes.
4. To turn shaker off, rotate the timer knob to 0 and the speed control to the extreme counterclockwise position. Depress the power switch to the off position.