

## **Centrifuges**

### **Sorvall RC5B (NSC 460, Kell 445)**

**Every time you use the centrifuge, you MUST sign the log book.**

**Balance all tubes.**

**Clean up any spills in the rotor and/or the centrifuge.**

Contact: Sonja Young (404) 413-5363; [sstovall@gsu.edu](mailto:sstovall@gsu.edu)

1. Turn power on. Front side panel on the right.
2. Door should already be open.
3. Put rotor in and tighten lid
4. Adjust speed knob.
5. Adjust time knob.
6. Adjust temperature knobs.
  - a. Blue needle – temperature setting
  - b. Red needle – max temperature setting
  - c. Black needle – actual temperature
7. Press start button.
8. When done, press door button on the front side panel, just above power button.
9. Leave door open.